


はくぶん  
HAKUBUN

# Air & Water Compression Experiment Kit - Model D

Made in Japan

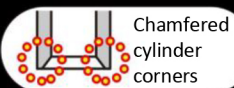

Chamfered  
cylinder  
corners

**Air compressor**

Easy to reassemble!  
You can easily change the position of the brim!

**Water compressor**

come with large  
syringe of 20ml

**Thick &  
Durable  
Center nozzle!**

Easy to close!

Comes with  
a bag clip

comes with  
a plastic bag

**Enrichment learning :  
fountain**

soft tube

Easy to fit!  
tube with a cap!

\*The bottle is  
not included in  
the set

**High screw  
thread  
prevent  
air escape!**


Sturdy, highly  
transparent cylinder!  
Made of special plastics!

**Stable  
design!**

Do not wipe off the lubricating oil "inside  
the cylinder" and "rubber of the pushrod".

## Attention!

- Before experimenting, please read the instructions carefully.
- Do not sit on top of the plastic bag filled with air.
- Do not put sponge balls or small parts in your mouth.
- Please do not wear a plastic bag on your head.
- Do not aim air or water compressors at people or windows.
- Please listen carefully to the teacher's instructions before using the tools.

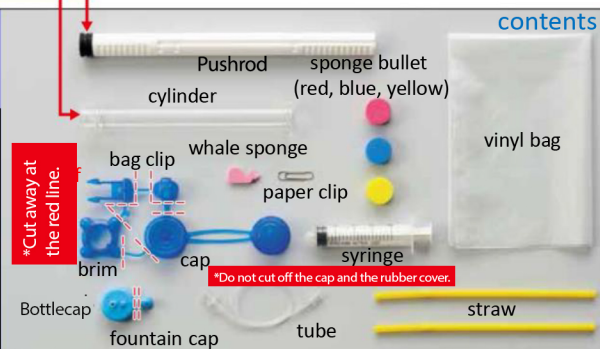

## Introduction: Let's collect some air

- Fill the plastic bag with air and check how it feels when holding it down.



Do not sit on top of the plastic bag. It is dangerous if torn,  
or the bag clip is blown off.

## How to trap air

Twist the top 15cm of the plastic bag three times, and insert the  
bag clip into the top 10cm of the bag until it clicks into place.

